

Enclosed Laboratory Filtration System BECO INTEGRA® LAB 60 P



Versatile, Sanitary Design Laboratory Depth Filtration System

BECO INTEGRA LAB 60 P from Eaton's Begerow Product Line is an enclosed single-sheet depth filtration system with a wide range of accessories for small volume filtration tasks in laboratories and research centers. The areas of application include small batches and upscaling in biotechnology and pharmaceutical, chemical, and cosmetics industries.

Special Features

- **Simple and clearly arranged design** with many connection, valve, and pressure measurement options.
- A wrap-around gasket ensures a **high level of safety** by preventing leakage to the surrounding atmosphere.
- Ergonomic design for **easy handling**. Effortless replacement of BECO® depth filter sheets; thorough and **easy cleaning**, no dead spaces.
- **Variable application possibilities** by using the base filter and sludge chambers of various size (with and without heating) for sheet and cake filtration.
- System components enable **parallel or step filtration** processes using one arrangement of several filters.

Other models and enhancements with additional components for complete systems are possible according to requirements.



Housing Material

Product contacting parts:	Stainless steel AISI 316L/AISI 316Ti
Others:	Stainless steel AISI 304
Gaskets, O-rings:	EPDM, Silicone, Viton, FEP coated

Connections

Filter inlet:	DN 8 ,Tri-Clamp, ISO 1127
Filter outlet:	DN 8 ,Tri-Clamp, ISO 1127

Technical Data


Max. permissible operating pressure:	87 psi (600 kPa/6 bar)
Max. permissible differential pressure:	58 psi (400 kPa/4 bar)
Max. permissible operating temperature:	284 °F (140 °C)
Depth filter size:	Ø 2.36", 0.08..0.16" thick (Ø 60 mm, 2...4 mm)

Pos.	Description	Article no.
10	BECO INTEGRA LAB 60 P (including 11.1, 11.2 (EPDM),12 + 13)	P7000302
12	Clamp BECO INTEGRA LAB 60 P	560630
13	Perforated plate BECO INTEGRA LAB 60 P	560604
40	Diaphragm pressure gauge 0 – 87 psi (0 – 600 kPa/0 – 6 bar), NG 63	P7400113
70	Adapter TC DN 8* - Hose clip**	P7500301
71	90° bend with 2 x TC DN 8*	P7500401
80	Cross piece 3 x TC DN 8*, 1 x TC DN 25***	P7500101
90	T-piece 2 x TC DN 8*, 1 x TC DN 25***	P7500201
100	Clamp DN 8	P7510101
120	Clamp DN 25	P7510001

Pos.	Description	8.45 oz (250 ml)	16.90 oz (500 ml)	33.8 oz (1,000 ml)
20	Filling space (without Pos. 11 + 12)	P7000602	P7000702	P7000802
30	Filling space with heating jacket (without Pos. 11 + 12)	P7000603	P7000703	P7000803

Pos.	Description	EPDM	Silicone	Viton	FEP coated
11.1	O-ring gasket BECO INTEGRA LAB 60 P for depth filters 0.08 – 0.14" (2 – 3.5 mm)	P7800305	P7800307	P7800306	P7800308
11.2	O-ring gasket BECO INTEGRA LAB 60 P for depth filters 0.14 – 0.18" (3.5 – 4.5 mm)	P7800309	P7800311	P7800310	P7800312
50	Diaphragm valve DN 8, 2 x TC DN 8*	P7400201	P7400202	P7400203	P7400204
60	Diaphragm valve DN 8, TC DN 8*/Hose clip**	P7400301	P7400302	P7400303	P7400304
110	Clamp gasket DN 8	P7510301	P7510302	P7510303	P7510304
130	Clamp gasket DN 25	P7510201	P7510202	P7510203	P7510204

* Tri-Clamp DN 8, ISO 1127 ** Hose clip DN 10 DIN 11854 *** Tri-Clamp DN 25, ISO 1127

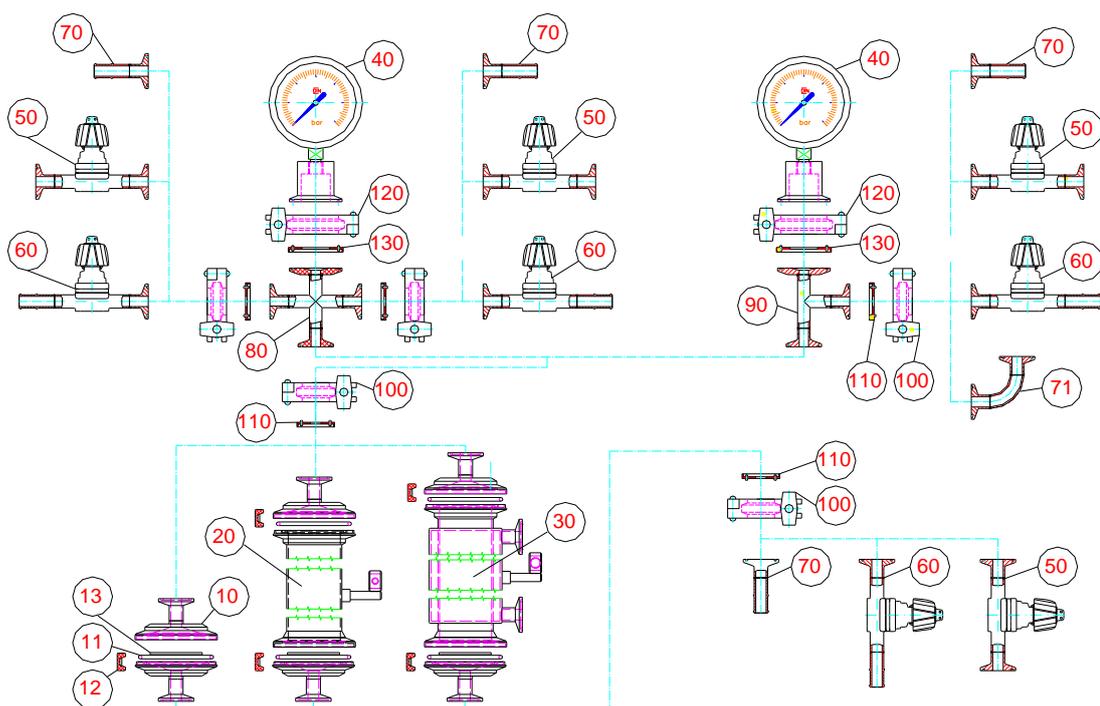

North America - HQ

44 Apple Street,
Tinton Falls, NJ 07724
Toll Free: 800 656-3344
(North America Only)
Voice: +1 732 212-4700

Europe/Africa/Middle East

Auf der Heide 2
53947 Nettersheim, Germany
Voice: +49 2486 809-0

Internormen Product Line

Friedensstraße 41
68804 Altlussheim, Germany
Voice: +49 6205 2094-0

Begerow Product Line

An den Nahewiesen 24
55450 Langenlonsheim, Germany
Voice: +49 6704 204-0

Brazil

Av. Julia Gaioli, 474 - Bonsucesso
07251-500 - Guarulhos
Brazil
Voice: +55 11 2465 8822

China

No. 7 Lane 280 Linhong Road,
Changning District,
Shanghai 200335, China
Voice: +86 21 5200 0422

Singapore

4 Loyang Lane #04-01/02
Singapore 508914
Voice: +65 6825 1668

For more information, please e-mail us at filtration@eaton.com or visit us online at eaton.com/filtration for a complete list of Eaton's filtration products.

Not all products in Eaton's Begerow Product Line are available in all regions. Please contact your local Eaton office to determine availability.

© 2012 Eaton Corporation. All Rights Reserved. All trademarks and registered trademarks are the property of their respective owners.

All information and recommendations appearing in this brochure concerning the use of products described herein are based on tests believed to be reliable. However, it is the user's responsibility to determine the suitability for his own use of such products. Since the actual use by others is beyond our control, no guarantee, expressed or implied, is made by Eaton as to the effects of such use or the results to be obtained. Eaton assumes no liability arising out of the use by others of such products. Nor is the information herein to be construed as absolutely complete, since additional information may be necessary or desirable when particular or exceptional conditions or circumstances exist or because of applicable laws or government regulations.