

BECO

Single Sheet Filter BECO INTEGRA[®] SOLO 300 – 1200

Universal Enclosed Depth Filtration System for a Diverse Range of Applications

The BECO INTEGRA SOLO single-sheet filter from Eaton's Begerow Product Line is ideal for difficult filtration tasks in the biotechnology, chemical, cosmetics, pharmaceutical, and food industries. The system separates and washes solids from suspensions after precipitation or crystallization processes, or for the processing of pre-coat residues.

The BECO INTEGRA SOLO mobile single-sheet filter system's filling space is designed as a pressure vessel with a level filter bottom that drops for cleaning, and a horizontal depth filter medium. The pressure vessel, with optional thermal/insulating jacket, closes with clamp screws and seals with an O-ring. BECO[®] depth filter sheet, filter cloth, or metal composite plate filtration is optional. For the manual discharge of the filter cake, the bottom part drops via a spindle or hydraulics.

Special Features

- **Multiple process steps** such as extraction, filtration, washing, dissolving, and discoloration in one device.
- Simple handling through ergonomic design. Easy replacement of the BECO depth filter sheet/filter cloth, simple and thorough cleaning
- **CIP** possible by using rotating spray nozzles that eliminate dirt residues in corners
- The design and construction of BECO INTEGRA SOLO single-sheet filter complies with the requirements of the GMP guidelines and the FDA regulations

Customization or additional components allow for system expansion or complete systems according to customer specifications.



Technical Data

Operating	-14.5/+87 psi
pressure:	(-100/+600 kPa, -1/+6 bar)
Operating	14 – 320 °F
temperature:	(-10/+160 °C)
Calculated temperature:	392 °F (200 °C)

Materials

Product contacting parts:	Stainless steel AISI 316L / AISI 316Ti Hastelloy, Enamel, fluor- polymeric coatings
Others:	Stainless steel AISI 304
Gaskets, O-rings:	EPDM, Silicone, Viton, FEP coated (FDA listed)
Filter material:	BECO depth filter sheet, filter cloth and/or metal compound plate
Connections:	Flange according to EN 1092-1 or tri-clamp, sterile flange, round thread etc.
Chassis:	Stainless steel AISI 304


Other possible features: agitator, quick closing device, tilting device, special designs possible







Туре	Depth filter sheets/depth filter media Ø in (mm)	Filter area ft² (m²)	Filling volume gal (I)	Solids capacity at a cake volume of 7.87 in (200 mm) gal (I)	Weight of filter Ib (kg)	А	Measur ii (m B		D
300	11.81	0.65	7.93	3.17	353	32.48	41.34	41.34	63.00
	(300)	(0.06)	(30)	(12)	(160)	(825)	(1,050)	(1,050)	(1,600)
450	17.72	1.72	17.17	8.45	551	38.39	41.34	47.24	55.12
	(450)	(0.16)	(65)	(32)	(250)	(975)	(1,050)	(1,200)	(1,400)
600	23.62	2.80	33.02	13.74	992	44.88	59.10	56.70	59.10
	(600)	(0.26)	(125)	(52)	(450)	(1,140)	(1,500)	(1,440)	(1,500)
800	31.50	5.17	84.54	23.36	2.425	54.33	53.15	70.08	78.74
	(800)	(0.48)	(320)	(96)	(1,100)	(1,380)	(1,350)	(1,780)	(2,000)
1000	39.37	8.18	166.43	40.15	3.748	63.78	66.93	82.68	80.71
	(1,000)	(0.76)	(630)	(152)	(1,700)	(1,620)	(1,700)	(2,100)	(2,050)
1200	47.24	11.73	290.59	57.59	5.512	73.23	72.83	92.52	83.46
	(1,200)	(1.09)	(1100)	(218)	(2,500)	(1,860)	(1,850)	(2,350)	(2,120)


Socket list	300 – 600	800	1000 – 1200	According to
Feed	DN 25	DN 40	DN 40	EN 1092-1
Filtrate	DN 25	DN 40	DN 40	EN 1092-1
Aeration/De-aeration			DN 25	EN 1092-1
Safety valve			DN 40	EN 1092-1
Safety fitting	DN 25	DN 25		EN 1092-1
Pressure gauge			DN 25	EN 1092-1
Stand by	DN 65	DN 65	DN 65	EN 1092-1

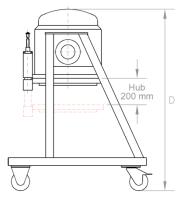

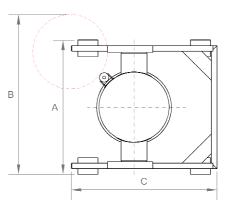

Single-sheet filter BECO INTEGRA SOLO 300 – 600

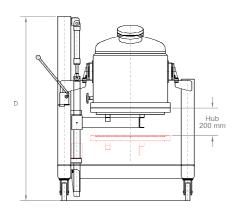

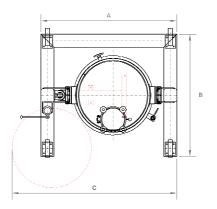

Single-sheet filter BECO INTEGRA SOLO 800 – 1200

North America - HQ

44 Apple Street, Tinton Falls, NJ 07724 Toll Free: 800 656-3344 (North America Only) Voice: +1 732 212-4700

Europe/Africa/Middle East

Auf der Heide 2 53947 Nettersheim, Germany Voice: +49 2486 809-0

Internormen Product Line Friedensstraße 41 68804 Altlussheim, Germany Voice: +49 6205 2094-0

Begerow Product Line An den Nahewiesen 24 55450 Langenlonsheim, Germany Voice: +49 6704 204-0

Brazil

Av. Julia Gaioli, 474 - Bonsucesso 07251-500 - Guarulhos Brazil Voice: +55 11 2465 8822

China

No. 7 Lane 280 Linhong Road, Changning District, Shanghai 200335, China Voice: +86 21 5200 0422

Singapore

4 Loyang Lane #04-01/02 Singapore 508914 Voice: +65 6825 1668 For more information, please e-mail us at filtration@eaton.com or visit us online at eaton.com/filtration for a complete list of Eaton's filtration products.

Not all products in Eaton's Begerow Product Line are available in all regions. Please contact your local Eaton office to determine availability.

© 2012 Eaton Corporation. All Rights Reserved. All trademarks and registered trademarks are the property of their respective owners.

All information and recommendations appearing in this brochure concerning the use of products described herein are based on tests believed to be reliable. However, it is the user's responsibility to determine the suitability for his own use of such products. Since the actual use by others is beyond our control, no guarantee, expressed or implied, is made by Eaton as to the effects of such use or the results to be obtained. Eaton assumes no liability arising out of the use by others of such products. Nor is the information herein to be construed as absolutely complete, since additional information may be necessary or desirable when particular or exceptional conditions or circumstances exist or because of applicable laws or government regulations.



EN 1 A 2.6.20.1 12-2012